


Assembly Instructions

Part Number: 21-808-0204

Product: PE 2000/3000
Torch/Spray Chamber
Adaptor Kit


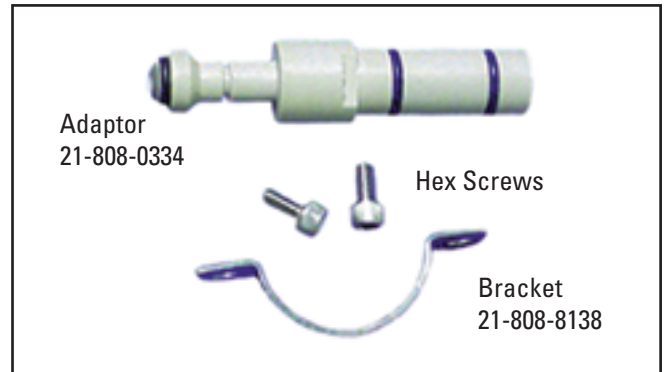


What You'll Need

A suitable injector, torch, spray chamber, nebulizer and ball joint clip.

Procedure

- 1) Insert injector into the adaptor.
- 2) Insert adaptor into torch
- 3) Insert assembled torch/adaptor/injector into mount block from the top of the mount block. The adaptor ball joint should just protrude through the bottom of the block exposing the groove/slot of the adaptor where the bracket should fit.
- 4) Position metal bracket in place and use metal screws to screw metal bracket to mount block.
- 5) Connect cup of spray chamber to torch adaptor ball joint, and hold in place with metal clip.

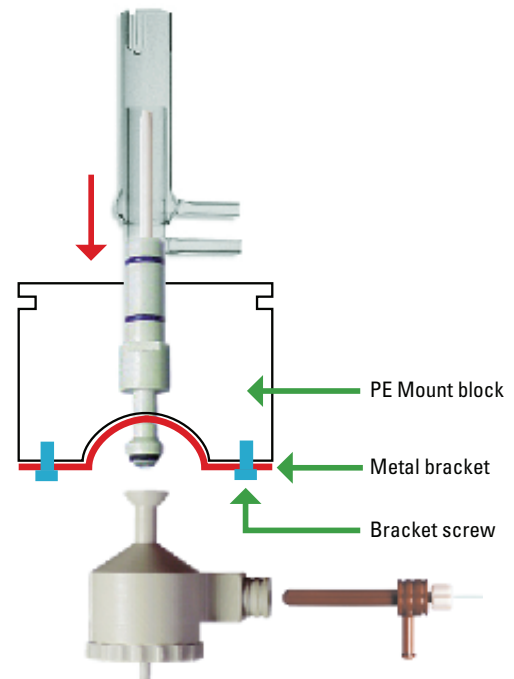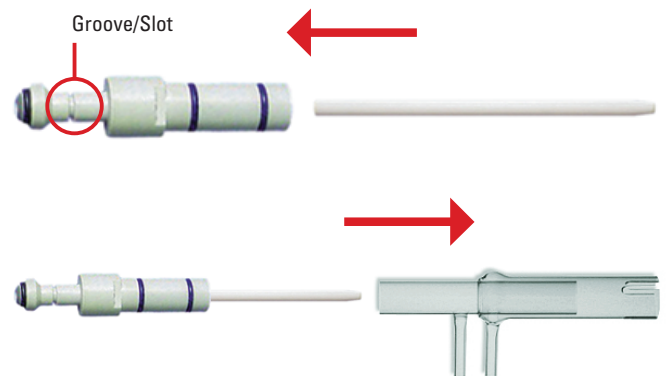

GLASS EXPANSION

Quality By Design

International

Glass Expansion
15 Batman Street
West Melbourne VIC 3003
Australia

Telephone: +61 3 9320 1111
Toll Free (Aust): 1800 777 638
Facsimile: +61 3 9320 1112
Email: enquiries@geicp.com

Americas

Glass Expansion
4 Barlows Landing Road
Unit 2A
Pocasset, MA 02559, USA

Telephone: 508 563 1800
Toll Free: 800 208 0097
Facsimile: 508 563 1802
Email: geusa@geicp.com

For your nearest Glass Expansion distributor, visit our website at www.geicp.com