


Assembly Instructions

Part Number: EL-2


Product: EzyLok Kit

1. P/N 6409-16: Tygon Tubing 34mm Piece
2. P/N 70-808-0735/2: 2 x EzyLok 6mm Hose Adaptor
3. P/N QSM-4: EzyLok Connector for 4mm Tubing
4. P/N SNP-1/6: 2 x Hose Clip 6mm

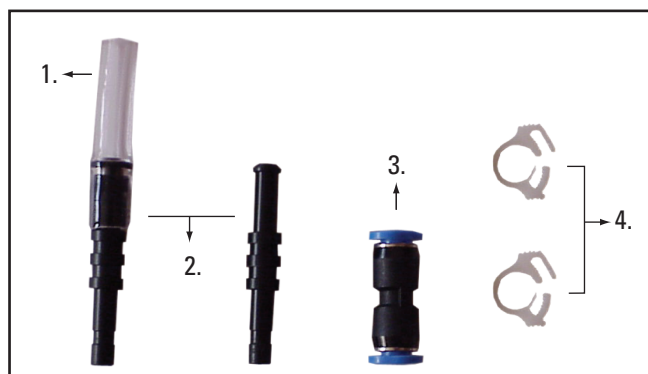


Note: To release or connect the EzyLok Connector, press down firmly on the blue button.

Using the EzyLok Kit with EzyLok Arm Nebulizers

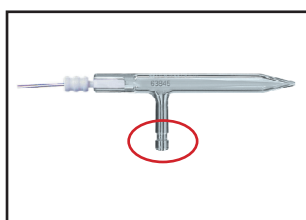


The EzyLok Arm

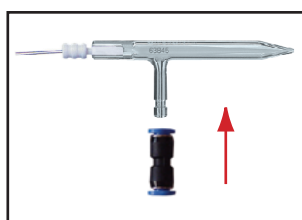


Step 1:

Connect the EzyLok Connector to the Nebulizer.

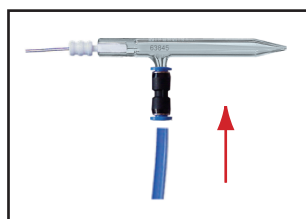

Step 2 (Option A)*:


If using Standard 4mm Tubing. Insert the tubing into the EzyLok Connector.

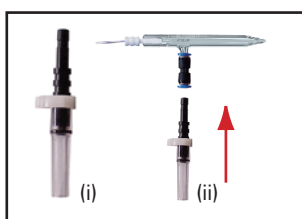


Step 2 (Option B)⁺:

If using other tubing. Secure tubing to Hose Adaptor with Hose Clip (i). The Clip should face the back of the nebulizer. Insert Hose Adaptor into Connector (ii).

Using the EzyLok Kit with Bulb Arm Nebulizers

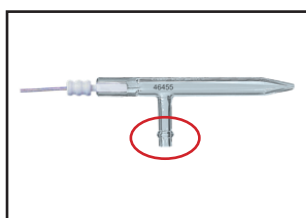


The Bulb Arm

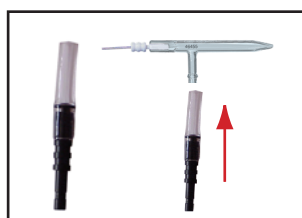

Step 1:

Push tube over bulb side arm until the tube meets the nebulizer body.

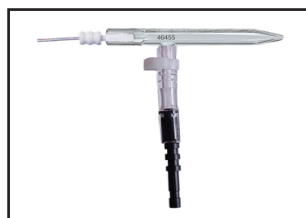

Step 2:

Secure tube with Hose Clip. Ensure Clip is between nebulizer glass body and bulb. Clip should face back of nebulizer.



Step 3:

Connect the EzyLok Connector to the Hose Adaptor.

Step 4: If using Standard 4mm tubing, continue by referring to *. If using other tubing, continue by referring to ⁺.

Important Note: Before applying gas pressure, apply a slight amount of pulling pressure between the Nebulizer and Hose to ensure they are securely locked.



GLASS EXPANSION

Quality By Design

International

Glass Expansion
15 Batman Street
West Melbourne VIC 3003
Australia

Telephone: +61 3 9320 1111
Toll Free (Aust): 1800 777 638
Facsimile: +61 3 9320 1112
Email: enquiries@geicp.com

Americas

Glass Expansion
4 Barlows Landing Road
Unit 2A
Pocasset, MA 02559, USA

Telephone: 508 563 1800
Toll Free: 800 208 0097
Facsimile: 508 563 1802
Email: geusa@geicp.com

For your nearest Glass Expansion distributor, visit our website at www.geicp.com