



Instructions for Inserting or Removing a Nebulizer


Hold the nebulizer by the barrel and slide it straight in or out.

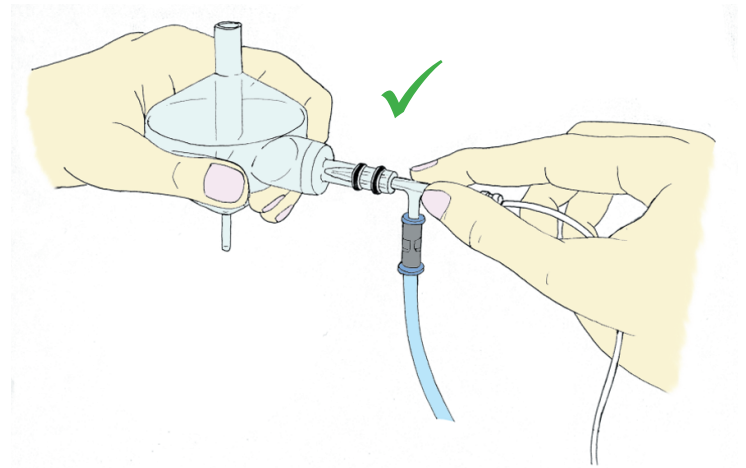

Do not hold the nebulizer by the side arm.

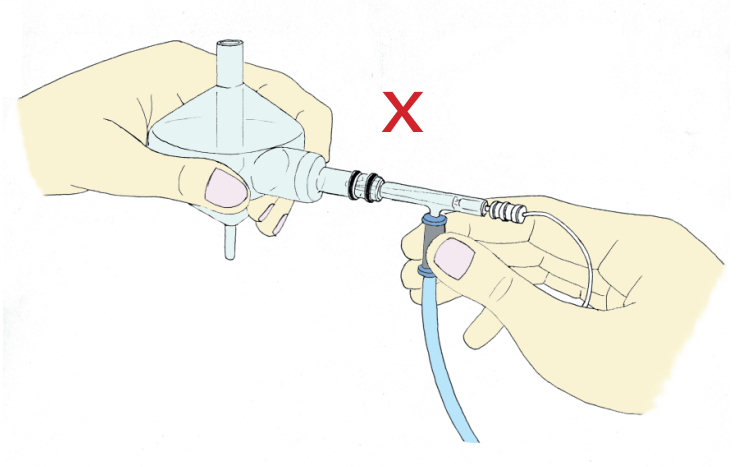

Do not apply any sideways force to the nebulizer. This may damage the nebulizer or the spray chamber.

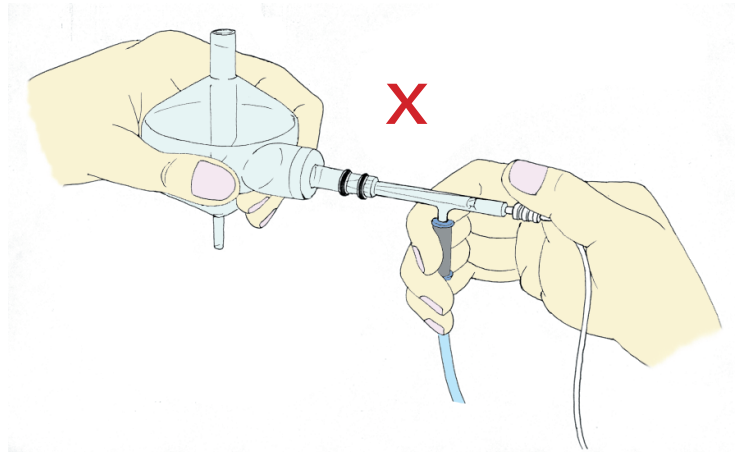

GLASS EXPANSION

Quality By Design

International

Glass Expansion
15 Batman Street
West Melbourne VIC 3003
Australia

Telephone: +61 3 9320 1111
Toll Free (Aust): 1800 777 638
Facsimile: +61 3 9320 1112
Email: enquiries@geicp.com

Americas

Glass Expansion
4 Barlows Landing Road
Unit 2A
Pocasset, MA 02559, USA

Telephone: 508 563 1800
Toll Free: 800 208 0097
Facsimile: 508 563 1802
Email: geusa@geicp.com

For your nearest Glass Expansion distributor, visit our website at www.geicp.com