

Running IsoMist-XR Software on Unsupported Operating Systems



Purpose

This guide assists customers facing difficulties with running the IsoMist-XR software on newer operating systems, such as Windows 11. If you're encountering issues with your IsoMist on a newer system, a common workaround is to try operating the software in compatibility mode. This document provides step-by-step instructions on how to do so.

Pre-Requisites

• The IsoMist-XR software must already be installed on your computer.

Instructions

These instructions are written for Windows 11, but the same steps apply to Windows 8.1 and Windows 10 operating systems at the time of writing.

1. Find the icon/executable for your IsoMist XR software. It's usually on your desktop or in the start menu.



2. Right-click the IsoMist XR icon and choose 'Properties'.



3. Click 'Change settings for all users' at the bottom of the window.



4. In the properties window, click the 'Compatibility' tab.

isolviist-Xit Pit	perces	
Security	Details	Previous Versions
General	Shortcut	Compatibility
If this program isn't try running the con	t working correctly on t npatibility troubleshoot	this version of Windows, er.
Run compatibil	ity troubleshooter	
How do I choose	compatibility settings m	anually?
Compatibility mod	de	
Run this prog	gram in compatibility mo	ode for:
Windows 8		~
Settings		
Reduced col	our mode	
8-bit (256) colou	ur 🗸	
🗌 Run in 640 x	480 screen resolution	
Disable full-se	creen optimisations	
Run this prog	gram as an administrato	pr
Register this	program for restart	
Use legacy d	lisplay ICC colour mana	agement
Change hi	igh DPI settings	
🔷 Change set	tings for all users	

5. Tick the box that says 'Run this program in compatibility mode for:'

Security	Details	Previous Versions	
General	Shortcut	Compatibility	
If this program isn't try running the cor Run compatibil How do I choose Compatibility mo S Run this prog	t working correctly on in npatibility troubleshoot ity troubleshooter compatibility settings n de gram in compatibility m	this version of Windows, er. nanually? ode for:	
Settings Reduced col 8-bit (256) color Run in 640 x Disable full-s	lour mode ur 480 screen resolution creen optimisations	1	
Run this prog	gram as an administrat	or	
Use legacy d	program for restart lisplay ICC colour man igh DPI settings	agement	
Change set	tings for all users		

- 6. The drop-down menu will now be enabled. Choose 'Windows 8'.
- 7. Click 'Apply' and 'OK'.
- 8. Open the IsoMist-XR software to check if it's working correctly.

For more information please visit **www.geicp.com** or contact one of our offices below.

Asia Pacific

6 Central Boulevard, Port Melbourne, VIC 3207, Australia Telephone: +61 3 9320 1111 Email: enquiries@geicp.com

Americas

31 Jonathan Bourne Drive, Unit 7 Pocasset, MA 02559, USA Telephone: 508 563 1800 Email: geusa@geicp.com

Europe

Friedenbachstrasse 9, 35781 Weilburg, Germany Telephone: +49 6471 3778517 Email: gegmbh@geicp.com