



GLASS EXPANSION

Quality By Design

Updating the Firmware on IsoMist XS




Purpose

This document provides a step-by-step guide on how to update the firmware on the IsoMist-XS unit using the IsoMist XS software.

Pre-Requisites

1. IsoMist-XS unit, with an active connection to the IsoMist XS software running on any computer.
 - The firmware update process is compatible with any operating system running the IsoMist software. This guide includes screenshots from both Windows and Linux software versions.

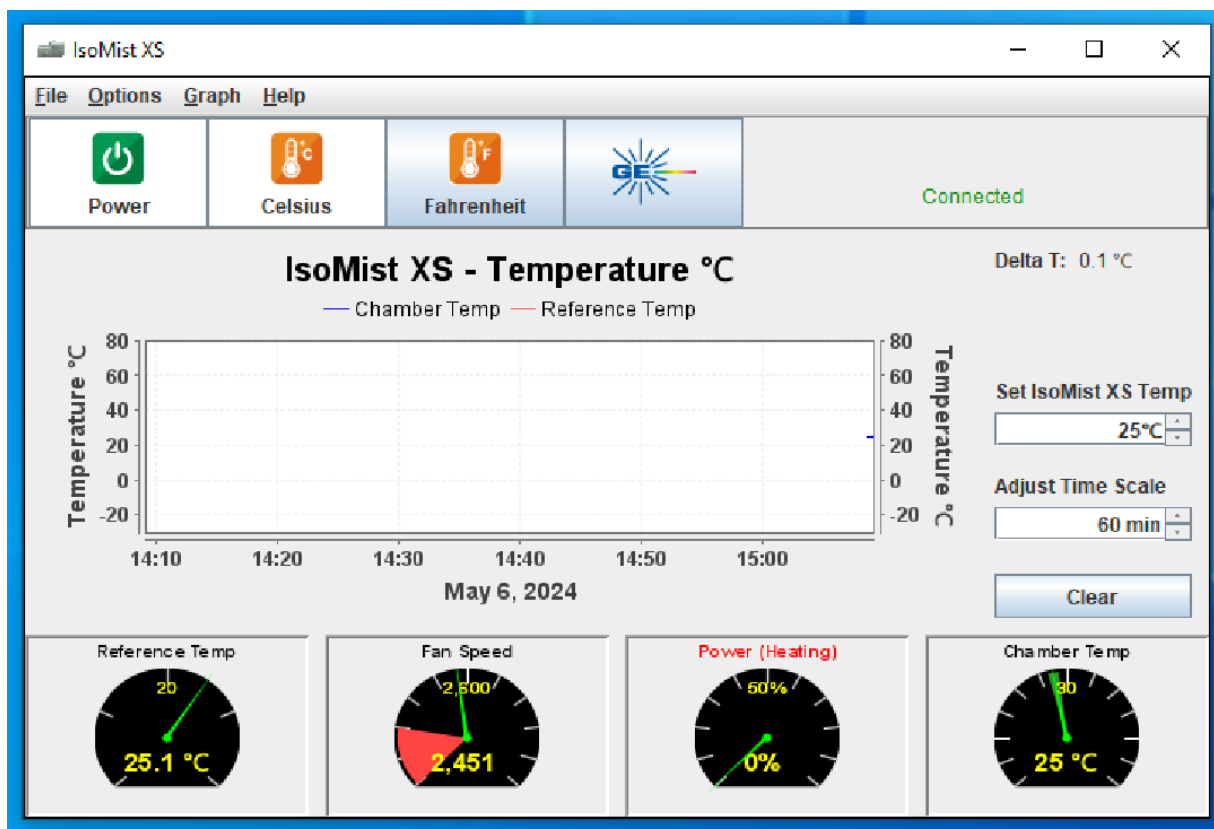


2. Copy of the IsoMist-XS firmware file.

- IsoMist firmware files use the .hex extension.
- The version number of the firmware is typically included in the file name.
 - Below is an example of the firmware file for v1.2:


Name	Date modified	Type	Size
isomistxs_1.2.hex	2/05/2024 12:33 PM	HEX File	2,027 KB

Instructions


1. Open the IsoMist XS software and navigate to **File > Upload Firmware**.

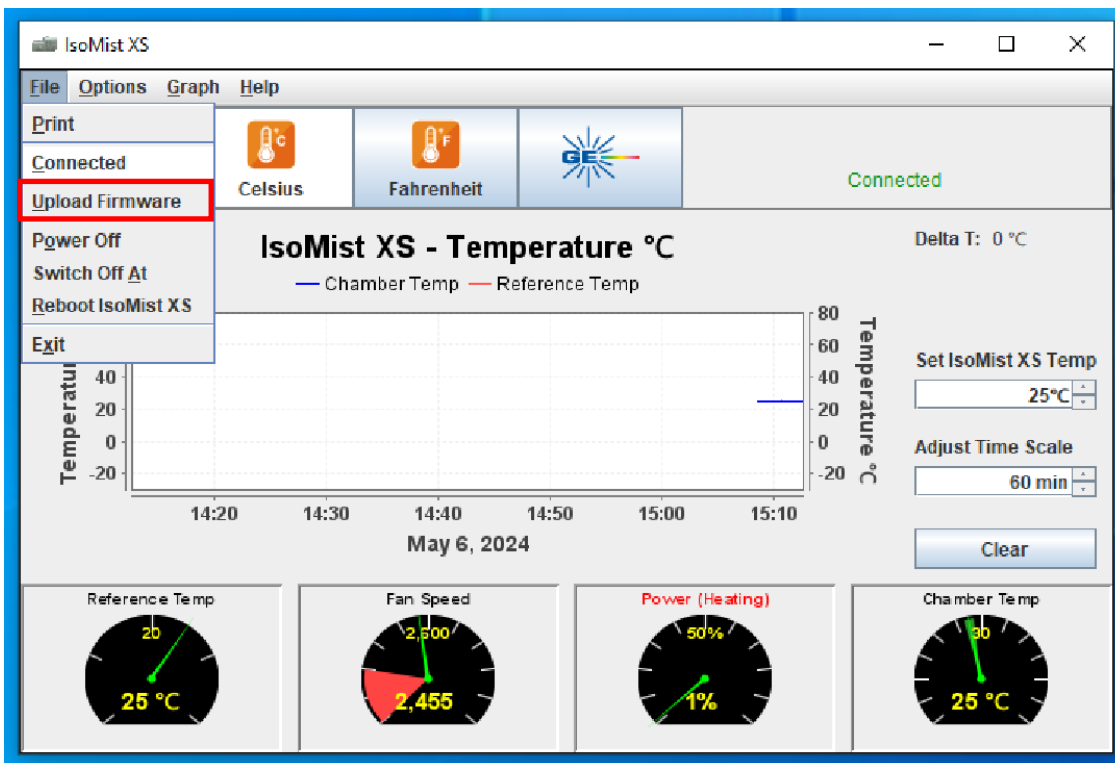


2. A pop-up will appear asking if you would like to apply a Firmware Update. Click **Yes**.

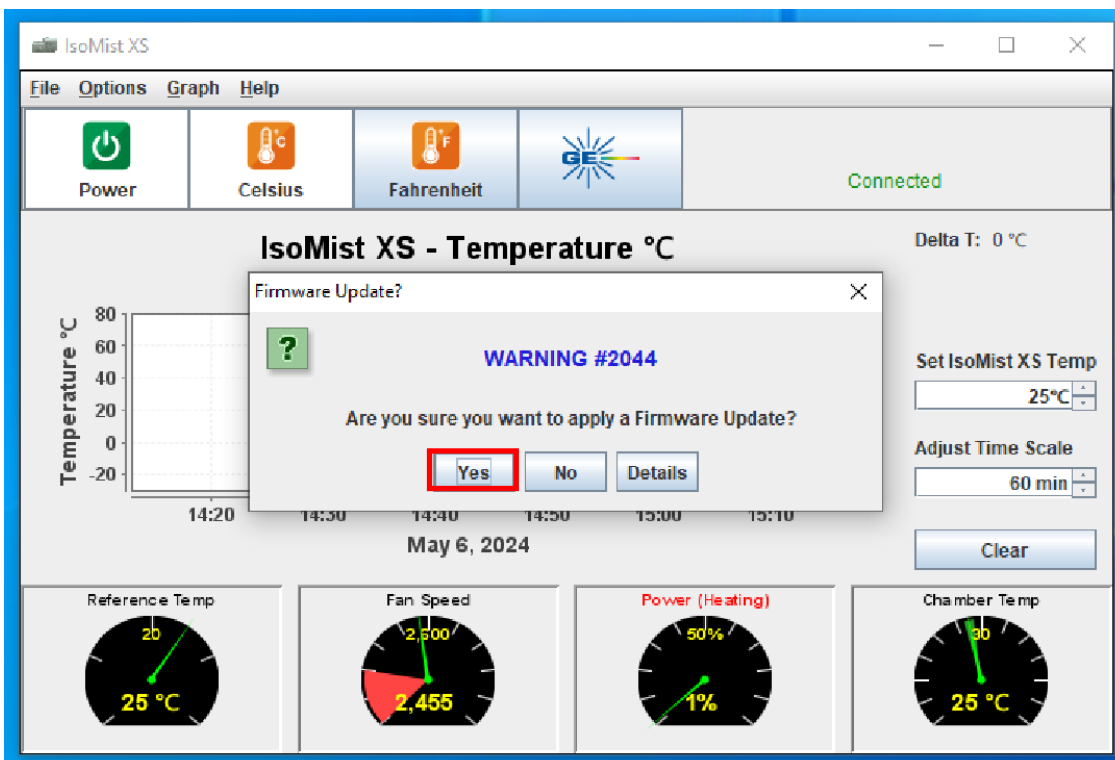

- The IsoMist will shut off briefly as it connects to the firmware. A notification will appear from the IsoMist during this process.

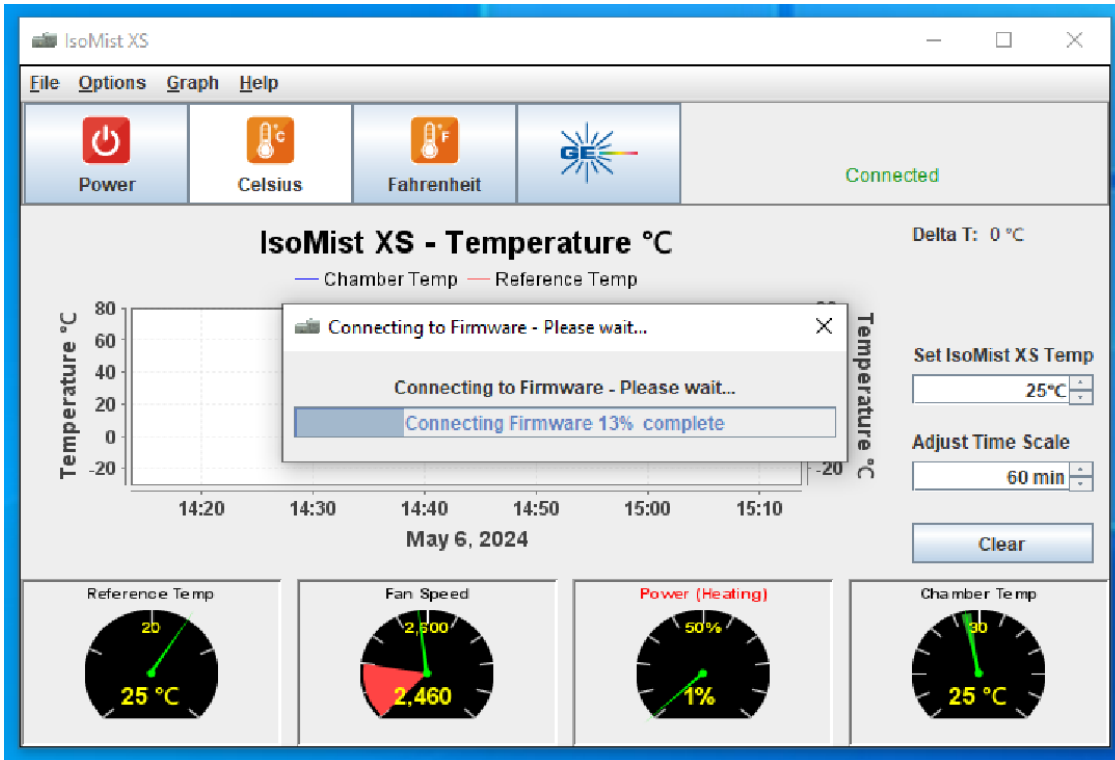


- Once the connection is established, a window will appear for you to select a firmware file. Choose the firmware file corresponding to the version you wish to upgrade to.

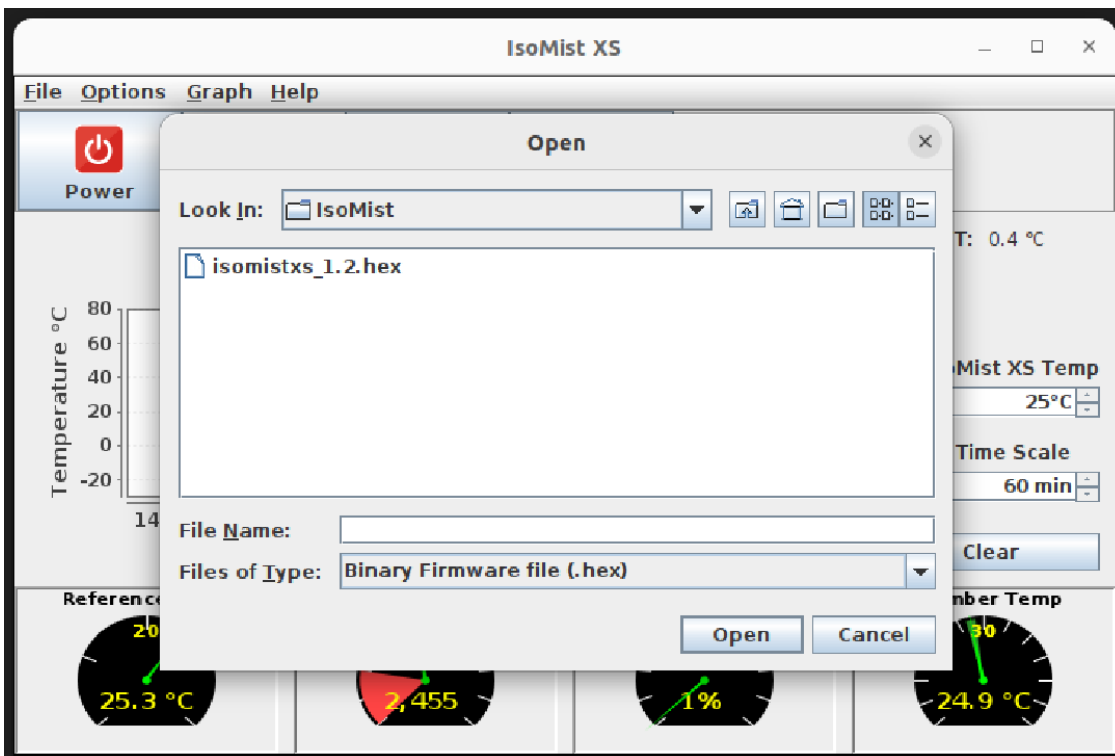



NOTE: If you are given the choice between multiple firmware files for a specific version, select the file with the **.hex** extension and ignore the **unified.hex** version.

- After selecting the firmware file, the application will begin the process of uploading the new firmware.

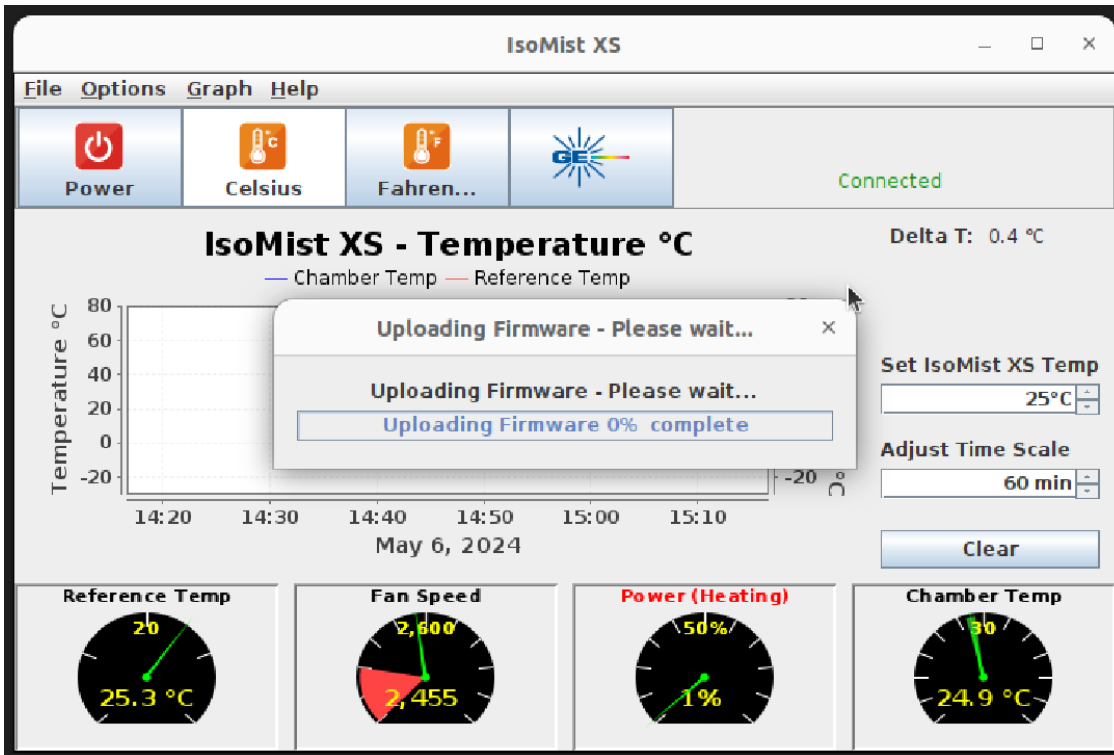

NOTE: The firmware update process may take 2-3 minutes. Do not turn off or disconnect the IsoMist during this step.

- When the update finishes, the IsoMist will automatically restart and re-establish a connection with the IsoMist software.
- To verify the firmware upgrade, navigate to **Help > About** in the IsoMist software and check if the firmware version matches the one you've updated to.



For more information please visit www.geicp.com or contact one of our offices below.

Asia Pacific

6 Central Boulevard, Port Melbourne, VIC
3207, Australia

Telephone: +61 3 9320 1111
Email: enquiries@geicp.com

Americas

31 Jonathan Bourne Drive, Unit 7 Pocasset,
MA 02559, USA

Telephone: 508 563 1800
Email: geusa@geicp.com

Europe

Friedenbachstrasse 9, 35781 Weilburg,
Germany

Telephone: +49 6471 3778517
Email: gegmbh@geicp.com