




# Replacing New Sample Tube & O'ring for PolyCon & OpalMist Nebulizers

1. Screw off sample tube thread cap by turning it anti-clockwise, and remove sample tube and o'ring from the nebulizer.

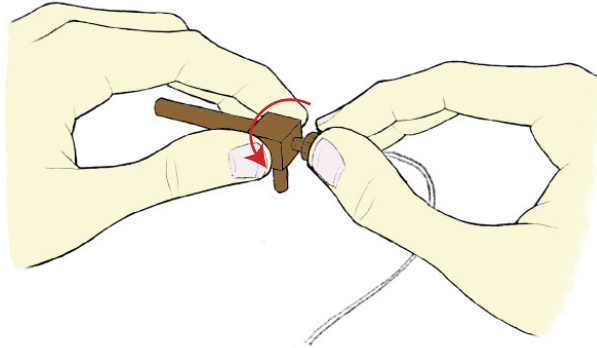

2. Cut a 45 degree angle on your **NEW** sample tube end by using a sharp blade, and insert it through the thread cap. Hold the end that you have just cut between your forefinger and your thumb. Make sure that the angle you have just cut is protruding no more than **2mm** between your thumb and forefinger. Place the new o'ring onto your forefinger, and insert the sample tube through the o'ring so that the o'ring sits about **25mm** away from the end that you have cut on an angle.

3. Before assembling the sample tube onto your nebulizer, **you must cut the end square** by using a sharp blade. Insert the sample tube into the nebulizer until it comes to a stop. Once you have come to a stop, grab hold of the thread cap and push it gently up towards the nebulizer. As you are pushing the thread cap it will make contact with the sample tube o'ring, keep pushing it forward until you reach the nebulizer. Once you have reached the nebulizer, start turning the thread cap clockwise until it comes to a stop.

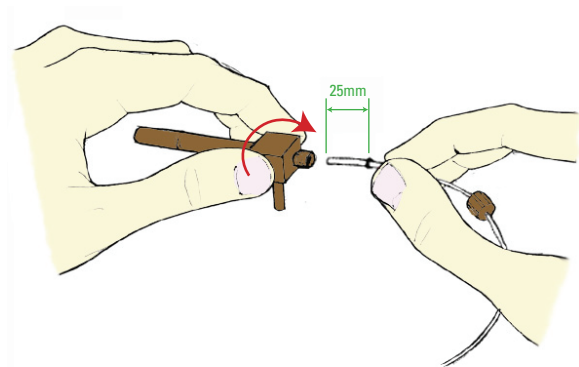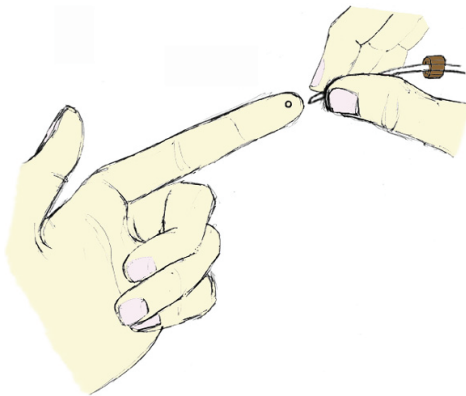

**Important Note:** Do not over tighten the thread cap, once you start to feel a bit of resistance with the thread cap, turn it another **1/4** of a turn and that should give you a reliable seal.



**Head Office (Australia)**  
15 Batman Street  
West Melbourne VIC 3003  
Australia  
A.B.N. 35 006 560 032

Telephone: +61 3 9320 1111  
Toll Free (Aust): 1800 777 638  
Facsimile: +61 3 9320 1112  
Email: [enquiries@geicp.com](mailto:enquiries@geicp.com)  
[www.geicp.com](http://www.geicp.com)

**USA Office**  
4 Barlows Landing Road  
Unit #2  
Pocasset, MA 02559  
USA

Telephone: 508 563 1800  
Toll Free: 800 208 0097  
Facsimile: 508 563 1802  
Email: [geusa@geicp.com](mailto:geusa@geicp.com)